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**ENGLISH**

Features ..... 05

Assembling and fitting instructions ..... 06

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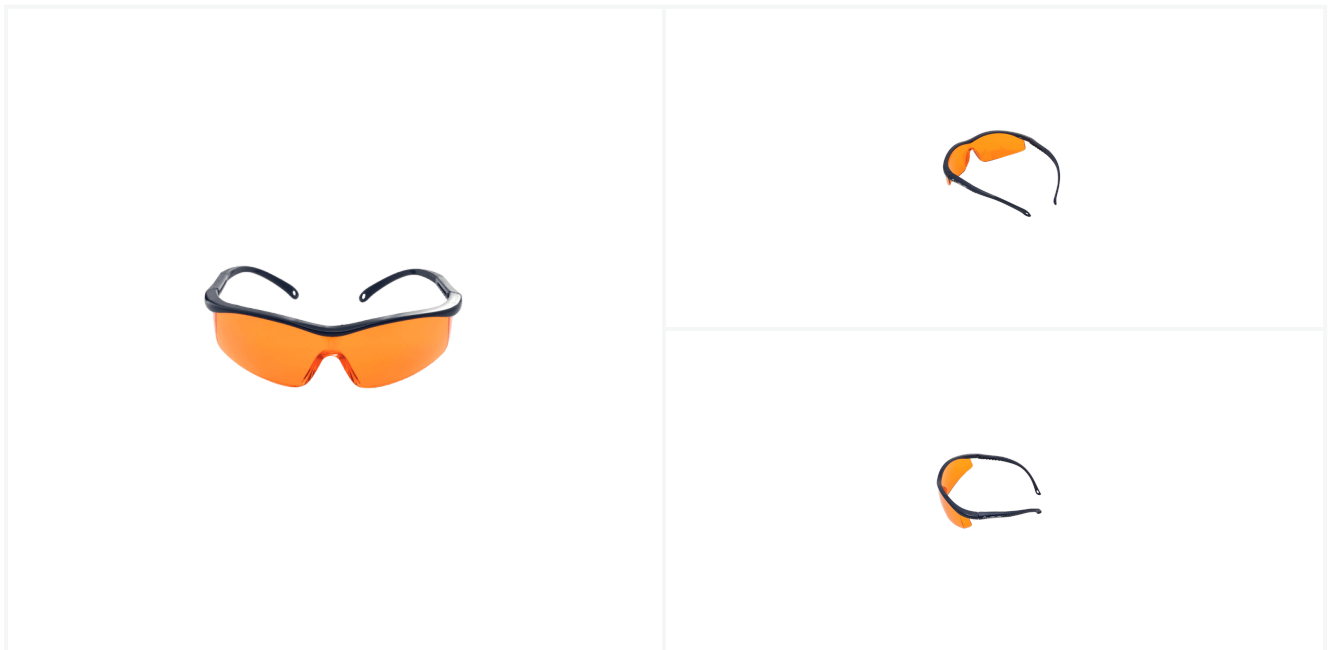
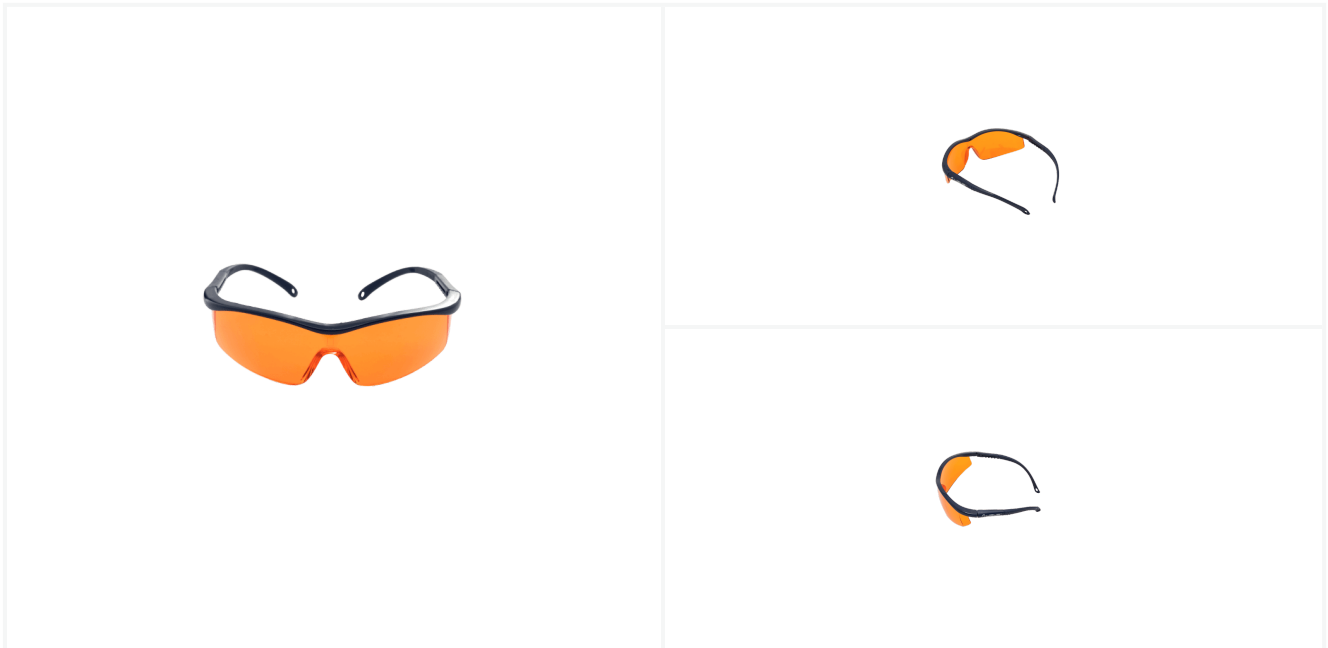
## FEATURES

- Designed to protect the eye against shock impact particles dust and sparks. Provides front and side protection.
- UV filter.
- Polycarbonate lens.
- Filters the violet and blue light present in the emission of UV lamps. Absorbs visible light up to 540nm. Applications: dentistry medicine laboratories UV curing of coatings and inks.
- Anti-fogging (AF): film resistant to condensation of moisture on the lens.

## FEATURES

- Stylized design of the ARGON model which makes it lighter and more ergonomic but with the same benefit.
- Flexible nylon frame and temples that provide a tight fit to the face.
- Telescopic arm adjustable to 4 length positions.
- Stainless steel screw.

## ASSEMBLING AND FITTING INSTRUCTIONS

### LIMITATIONS AND WARNINGS

1. It is made of materials that are not expected to cause adverse health effects; however contact with the user's skin may cause allergies in highly sensitive individuals. If the protector is damaged it should be discarded and replaced. This goggle provides the protection indicated in the technical specifications and according to its marking. Eye protectors marked with the letter "S" should not be used when there is a risk of impact from hard sharp or pointed flying particles. Operating temperature: -5°C to +40°C.

### USE

1. The life of the goggle once removed from its original packaging will depend on the frequency and conditions of use. For daily use it is recommended to change it every 6 months but under extreme conditions the protector may deteriorate faster; in either case discard the product and replace it with a new one. Do not attempt to repair it. No replacement parts are provided for this product. This protector should not be used over normal prescription eyeglasses. If the goggle is damaged during use immediately leave the hazardous area and replace it with a new one.

## CLEANING INSTRUCTIONS

1. Recommended after each use with warm running water and neutral soap or a cleaning solution specified for lenses. Do not use degreasing or chlorinated fluids or solvents.

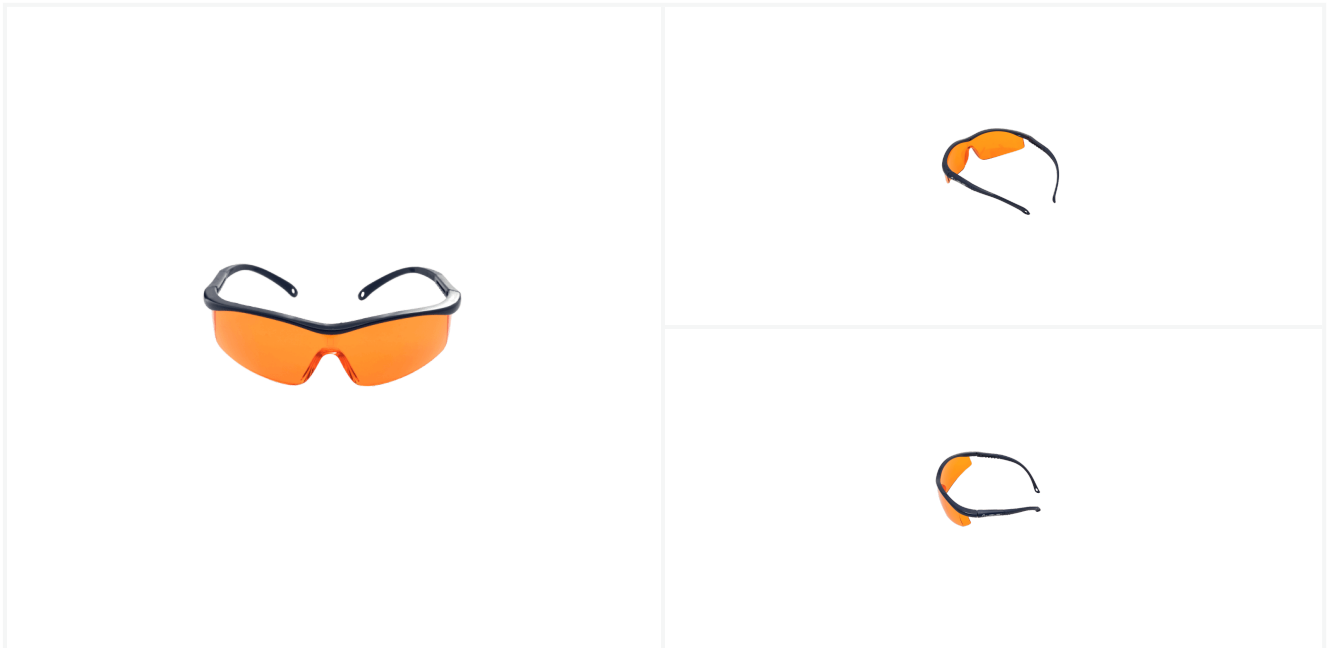
## **CLEANING INSTRUCTIONS**

### **STORAGE**

1. Keep in original packaging in dry clean place away from direct sunlight high temperature sources vapors solvents or chemicals. Temperature: -5°C to +50°C / Ambient relative humidity < 80%.

## **SPECIFICATIONS MET**

- EN 166
- ANSI Z87.1.
- You can know in detail the certified codes under the different international standards by consulting them on our website [www.libus.com](http://www.libus.com) and downloading the corresponding documentation accompanying the licenses.

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